


Starting to Work on the QuantStudio1 and QuantStudio3

For new users, please create a new user here- [BookitLab](#).

For booking time slots on the instruments, you will need to schedule training with the ATGC. For the QuantStudio1 [QS1] in the Emerson building, please contact [Nitsan](#). For the QuantStudio3 [QS3] in the Rappaport building, please contact [Liza](#).

Making a reservation

To reserve the instrument, go to INSTRUMENTS-calendar on the left menu. In the search box type 'Real Time PCR QS1 Emerson' or 'Real Time PCR QuantStudio3' to open the instrument's calendar.

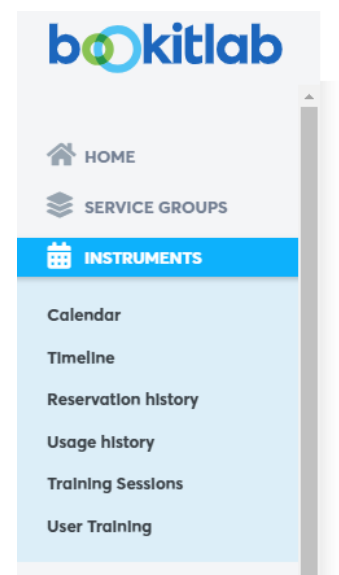

Start use

To start your reservation on QS1 or QS3, please login into your BookitLab account on the instrument's computer (the BookitLab software is running in the background).

BOOKITLAB
CORE FACILITIES MANAGEMENT SOLUTION

P4b003


User Name:

Password:

Login

⌵ Details:

After login, activate your booking-

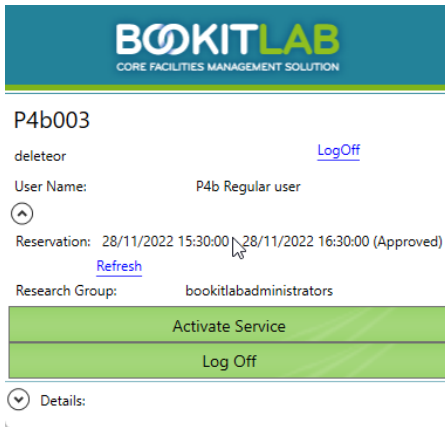

BOOKITLAB
CORE FACILITIES MANAGEMENT SOLUTION

P4b003

deleteor [Log Off](#)

User Name: P4b Regular user

Reservation: 28/11/2022 15:30:00 - 28/11/2022 16:30:00 (Approved)
[Refresh](#)

Research Group: bookitlabadministrators

Activate Service

Log Off

Details:

Your booking for the QuantStudio system will automatically end when your reservation time concludes. If you need to end your session earlier, log in to the BookitLab software and select the "Log Off" option. **Be aware that the QuantStudio software will close either when your reserved time ends or immediately after you log off. To avoid disruptions, ensure that you reserve enough time in advance to complete your work.**

Please note: Data files are stored in user-specific folders on the computer and can be accessed directly without requiring the QuantStudio software.

Should you require guidance or assistance, please do not hesitate to reach out.

ATGC team.

Emerson lab- 073-3781387 ; Rapport lab- 073-3785221.